

SAMPLE SUBMISSION GUIDE

Vector for Subcloning

Sending samples according to the requirements below helps us to do our job better and provides you with accurate results.

Sample Requirements

- Please send us all samples at a concentration of 200ng/μl with a 260/280 ratio of approximately 1.8.
- The total sample volume should be minimum 10μl (min. amount: 10μg)

If many subclonings have been ordered into the same vector, please send a correspondingly larger amount of vector (in one tube):

| | |
|--|----------|
| 2-5 Subclonings | 10-15 μg |
| 6-10 Subclonings | 15-20 μg |
| additional 5 subclonings please add 5 μg DNA | |

- Please do not perform a restriction digestion of your vector beforehand
- The sample should be sent in doubled distilled water or TE buffer
- Please send your samples in 1.5ml tubes
- Labelling of the tubes:
 - Sample Name as indicated in the request form
 - Concentration of the sample
 - Volume of the sample

Shipping

- Please mention the Order ID
- Please use a padded envelope when sending the vector to ensure that it arrives safely
- Ship the samples liquid at ambient temperature
- Shipping address:

Gene Synthesis - Lab
Eurofins Genomics
c/o Sarah Ebner
Anzinger Str. 7a
D-85560 Ebersberg

Contact

Please do not hesitate to contact us in case of any questions!

Email: genes-eu@eurofins.com; Phone: +49 7531 816 068

